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Cooling water

Certain requirements are placed on the cooling water with regard to purity. Depending on the cooling water contamination, a suitable method of purification and/or treatment of the water must be employed. The condenser and the complete cooling water circuit can become blocked, damaged and leaky due to unsuitable cooling water. Extensive consequential damage may arise on the whole cooling circuit. The cooling water quality depends on local conditions. If a fault or damage occurs due to unsuitable water quality, it is not covered by our guarantee.

Important: Danger of corrosion of the cooling water circuit due to water of unsuitable quality.

- Free chlorine (e.g. from disinfectants) and water containing chlorine lead to pitting in the cooling water circuit.
- Distilled, deionized or demineralized water is unsuitable due to its corrosive properties and leads to corrosion in the cooling water circuit.
- Seawater is unsuitable due to its corrosive properties and leads to corrosion in the cooling water circuit.
- Water containing iron or iron particles leads to rust formation in the cooling water circuit.
- Due to the high lime content hard water is not suitable for cooling and leads to calcification in the cooling water circuit.
- Cooling water with suspended matter is not suitable.
- Untreated and unpurified river or cooling tower water is not suitable due to its microbiological content (bacteria), which can become deposited in the cooling water circuit.
- Putrid water is not suitable.

Suitable cooling water quality

pH – value	7.5 – 9.0
Sulfates [SO ₄ ²⁻]	< 70 mg/L
Hydrocarbonates [HCO ₃ ⁻]/ sulfates [SO ₄ ²⁻]	> 1.0
Total hardness	4.0 – 8.5 °dH
Hydrocarbonates [HCO ₃ ⁻]	70 – 300 mg/L
Conductivity	10 - 500 µs/cm
Chlorides (Cl ⁻)	< 50 mg/L
Sulfites [SO ₃ ²⁻]	< 1 mg/L
Free chlorine gas (Cl ₂)	< 1 mg/L
Nitrates (NO ₃ ⁻)	< 100 mg/L
Ammonia (NH ₃)	< 2 mg/L
Iron (Fe), dissolved	< 0.2 mg/L
Manganese (Mn), dissolved	< 0.1 mg/L
Aluminum (Al), dissolved	< 0.2 mg/L
Free aggressive carbonic acid (CO ₂)	< 5 mg/L
Hydrogen sulfide (H ₂ S)	< 0.05 mg/L
Algae growth	Not permissible
Suspended matter	Not permissible

Risk to the environment due to oil contamination of the cooling water circuit

With a leaky condenser there is the danger that refrigerating machine oil from the coolant circuit of the cooling thermostat can pass into the cooling water.

Follow all the legal requirements and the regulations of the water supply utility which apply at the point of use.

Water pollution due to leakage

To avoid pollution due to a leak in the cooling water system it is recommended that a leakage-water detector with a water cut-off is installed.

Servicing intervals

Cleaning the dirt trap


At regular intervals of one month or longer, the dirt trap must be cleaned, depending on the degree of soiling.

Decalcifying the water cooling circuit

At regular intervals of 3 months or longer, the water-cooled condenser must be decalcified or cleaned. This depends on the hardness of the cooling water and the degree of soiling.

Required equipment:

- Two containers of approx. 10 to 20 liters volume.
- Use a suitable pump (drum pump) or a hose with funnel. Place the funnel as high as possible so that the device can fill quickly.
- Fit connecting hoses between container, pump, cooling water inlet and between cooling water outlet and back to container.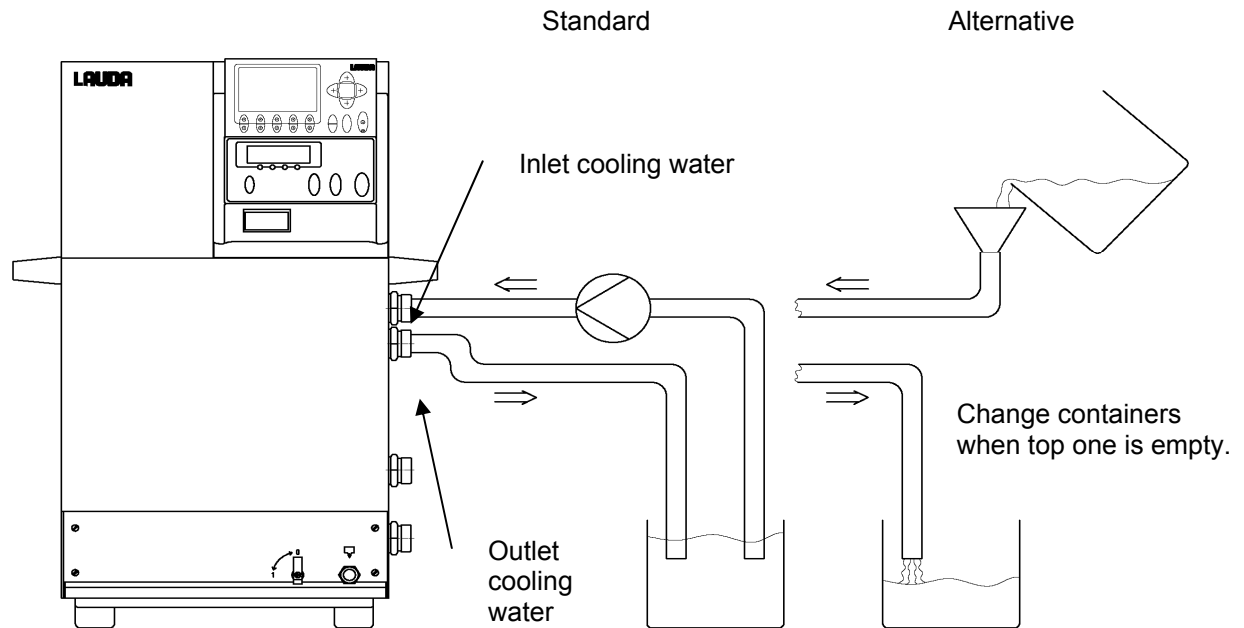

Via the water inlet hose, fill the device with decalcifier (pump or hose). To do this, set the setpoint to 10 °C. After starting the compressor, the water circuit can be filled. Circulate the decalcifier with the pump resp. continue to top up the decalcifier as necessary. Allow the decalcifier to take effect (refer to table below). Drain the device. Reconnect the device to the water supply and thoroughly flush out (refer to table below). While liquid is being pumped around the cooling water circuit, operate the device as described above at 10 °C.

Acting time	Continue the pump stage until most of the foamy reaction, usually at the start, has decayed. Generally, this is achieved after about 15 to 30 minutes.
Decalcifier	Water with LAUDA Decalcifier LZB 126 (5 kg). It is essential to follow the safety instructions and the handling instructions at the packing when handling the chemicals.
Flushing:	Allow at least 30 liters of water to flow through.

For Integral XT refer to special information in the Operating Instructions.