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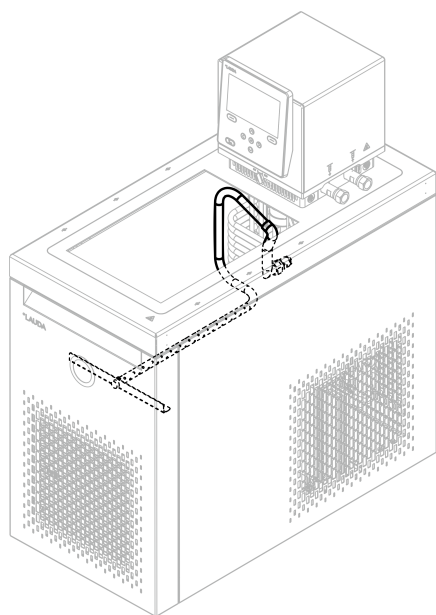
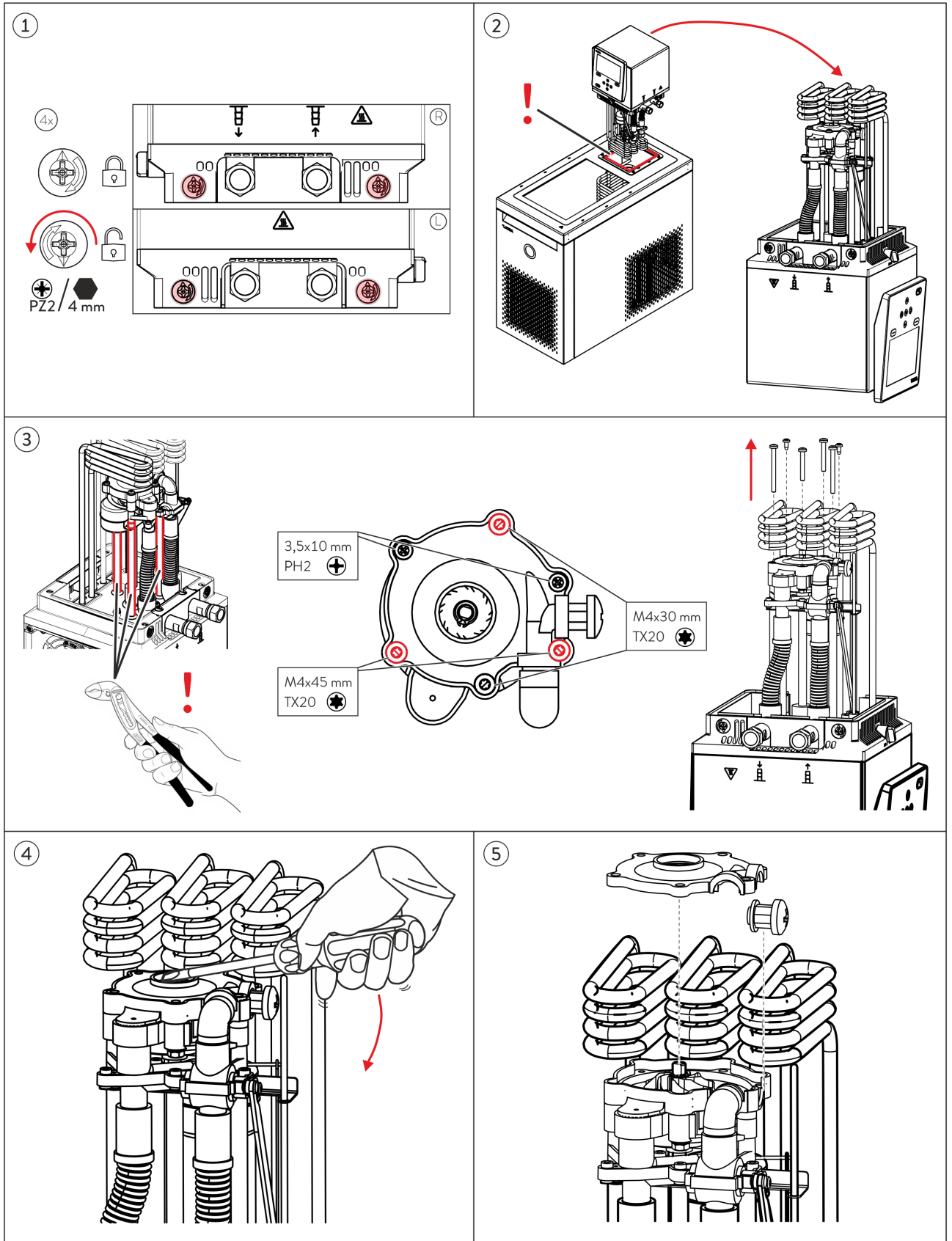
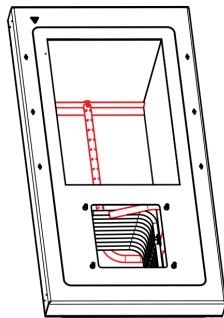
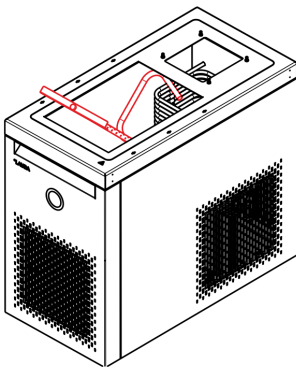
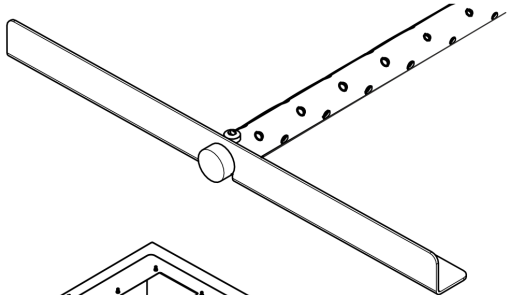
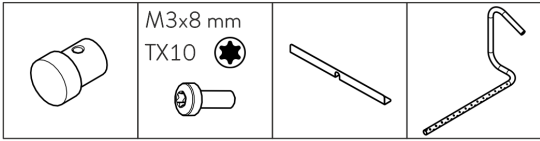
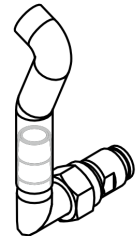
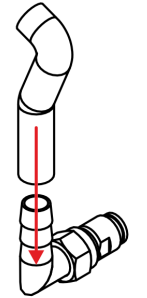
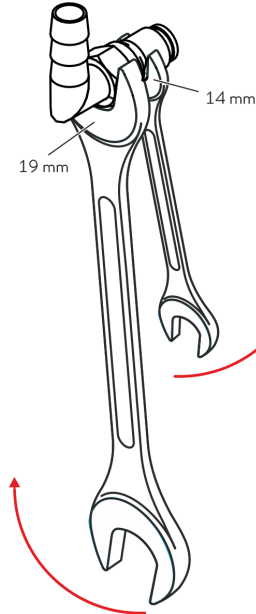
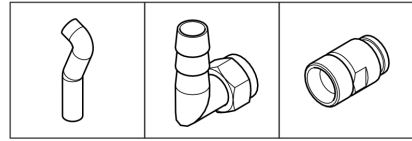
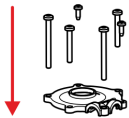
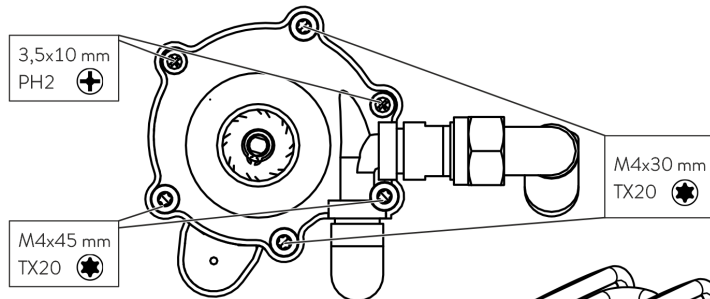
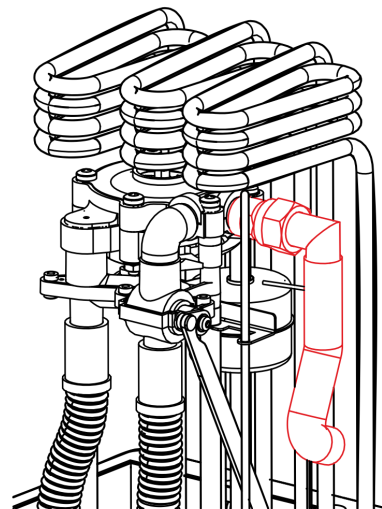
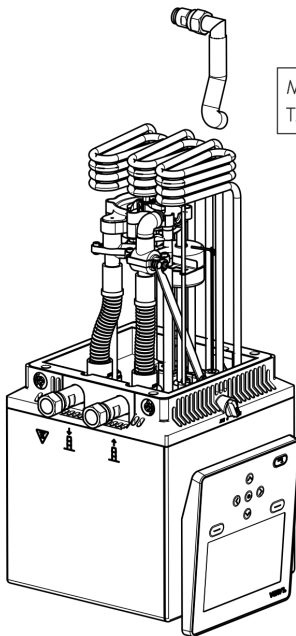
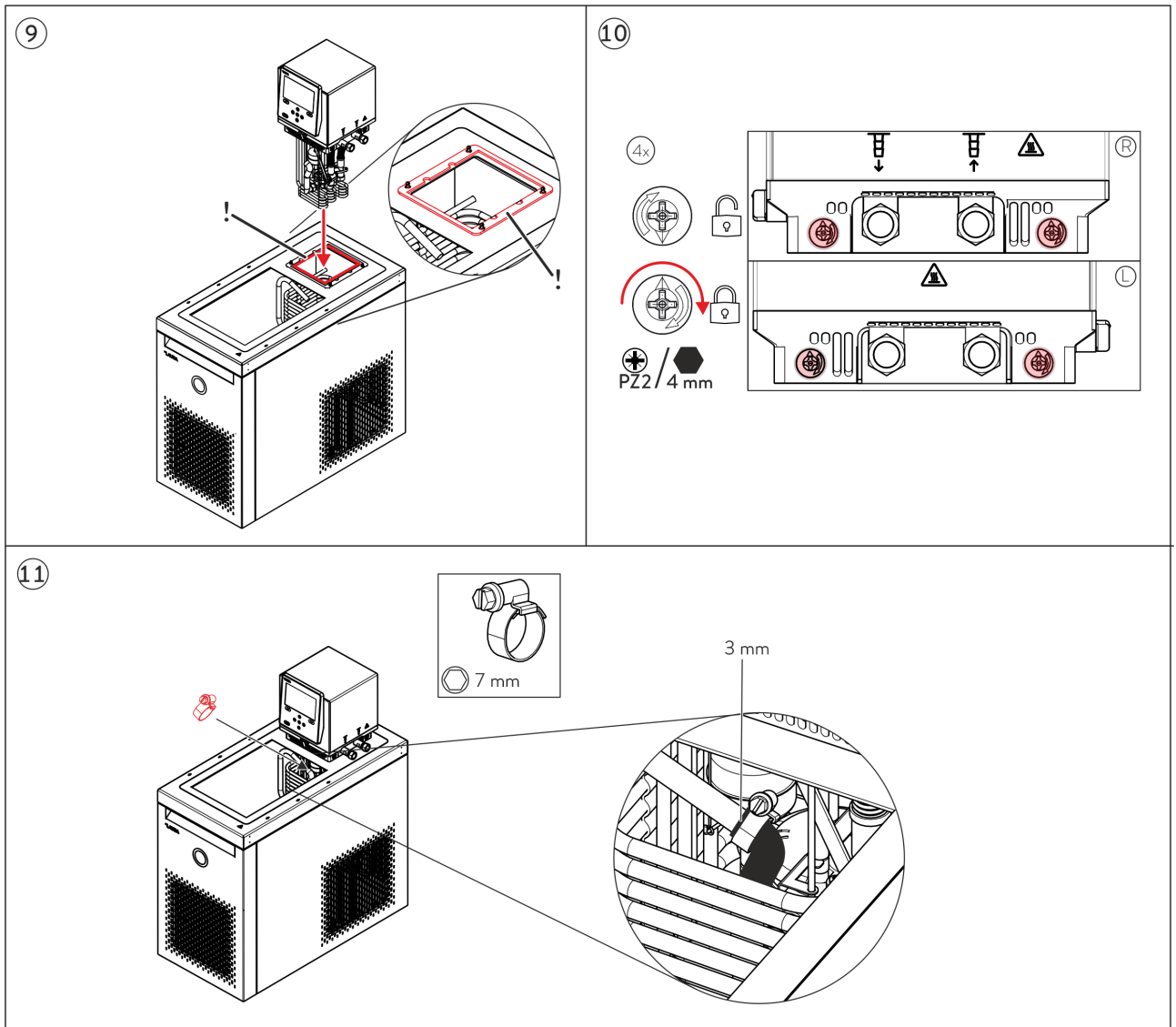
M4x45 mm
TX20

1 Introduction

These instructions describe how to mount a jet pipe on a LAUDA Universa MAX bath thermostat and apply to the following accessory sets. Read the instructions carefully before starting work.

Order number	Description
A001785	Jet pipe suitable for U 4230 M
A001786	Jet pipe suitable for U 1645 M, U 16 M
A001827	Jet pipe suitable for U 12 M, U 1245 M
A001828	Jet pipe suitable for U 20 M, U 2040 M

Intended use

This accessory set may only be used under the conditions specified in the operating instructions for the device. Any other mode of operation is considered improper use. It is the operator's responsibility to ensure that the device is used properly.

Applicable documents

- Operating manual Universa MAX and Universa PRO (Document number Q4DT-E_13-020)

For all work on the device that is necessary due to the mounting of the jet pipe and depending on the condition of the device, the instructions, warnings, and safety instructions in the operating manual must be observed.

Warranty conditions and disclaimer

LAUDA grants a standard warranty of one year. The components have been manufactured in accordance with high quality and safety requirements. LAUDA DR. R. WOBSE GMBH & CO. KG warranty and liability does not cover personal injury or property damage that can be attributed to one or more of the following causes, for example:

- Failure to follow these mounting instructions
- Improper mounting, commissioning, maintenance, and operation
- Improperly performed repairs and transport
- Unauthorized structural modifications to the device and accessories
- Any damage caused by continued use despite an obvious defect.
- No use of original spare parts and original accessories.
- Improper use of the device and accessories.

Personnel qualification

Specialized personnel

Certain activities on the device must be performed by specialized personnel. Specialized personnel are people whose professional education, knowledge, and experience as well as knowledge of relevant standards qualify them to assess the function and risks associated with the device and its use.

Personal protective equipment

Protective work clothing

Protective clothing is required for the work. This must meet the legal requirements for personal protective equipment. The protective clothing should be long-sleeved. Do not wear rings, chains or other jewelry.

Safety gloves

The safety gloves are used to protect against injury when removing the outer housing panels and when removing components. The safety gloves must be chemical-resistant for work involving possible contact with temperature control fluid.

Safety shoes

Safety shoes are used to protect against heavy falling parts and slipping on slippery surfaces. They are also used to protect the feet when removing the outer housing panels..

Structure of warnings

Warning signs	Type of danger
	Warning – hot surface.
	Warning – danger zone.
Signal word	Meaning
WARNING!	This combination of symbol and signal word indicates a potentially dangerous situation that can result in death or serious injury if it is not avoided.
CAUTION!	This combination of symbol and signal word indicates a possible dangerous situation that can result in minor injury if it is not avoided.

2 Safety instructions

WARNING!
Contact with live parts

Electric shock

- Disconnect the device from the power supply before performing any mounting work.

WARNING!
Danger of the device rolling away or overturning

Impact, crushing

- Do not tilt the device!
- Position the device on an even, non-slip surface with a sufficient load carrying capacity!
- Actuate the caster brake when setting up the device!
- Do not place heavy parts on the device!

WARNING!
Incorrect handling

Burning, fire, device damage

- Do not damage the cooling circuit.
- Do not use aggressive cleaners to clean the pump and control unit.
- Do not use chlorine-based cleaners for the bath vessel and evaporator.
- Do not use sharp or pointed objects to clean the evaporator.

CAUTION!
Risk of hand and foot injuries!

When installing the jet pipe, there is a risk of hand injuries from sharp edges and crushing, as well as foot injuries when removing and attaching the pump and control unit.

- Geeignete Schutzhandschuhe, Arbeitsschutzkleidung und Sicherheitsschuhe tragen.
- Wear suitable protective gloves, protective work clothing, and safety shoes.
- Ensure that the work area is clean, well lit, and free of obstacles.
- Always position yourself and work in such a way that your hands and feet are not below the parts to be removed.

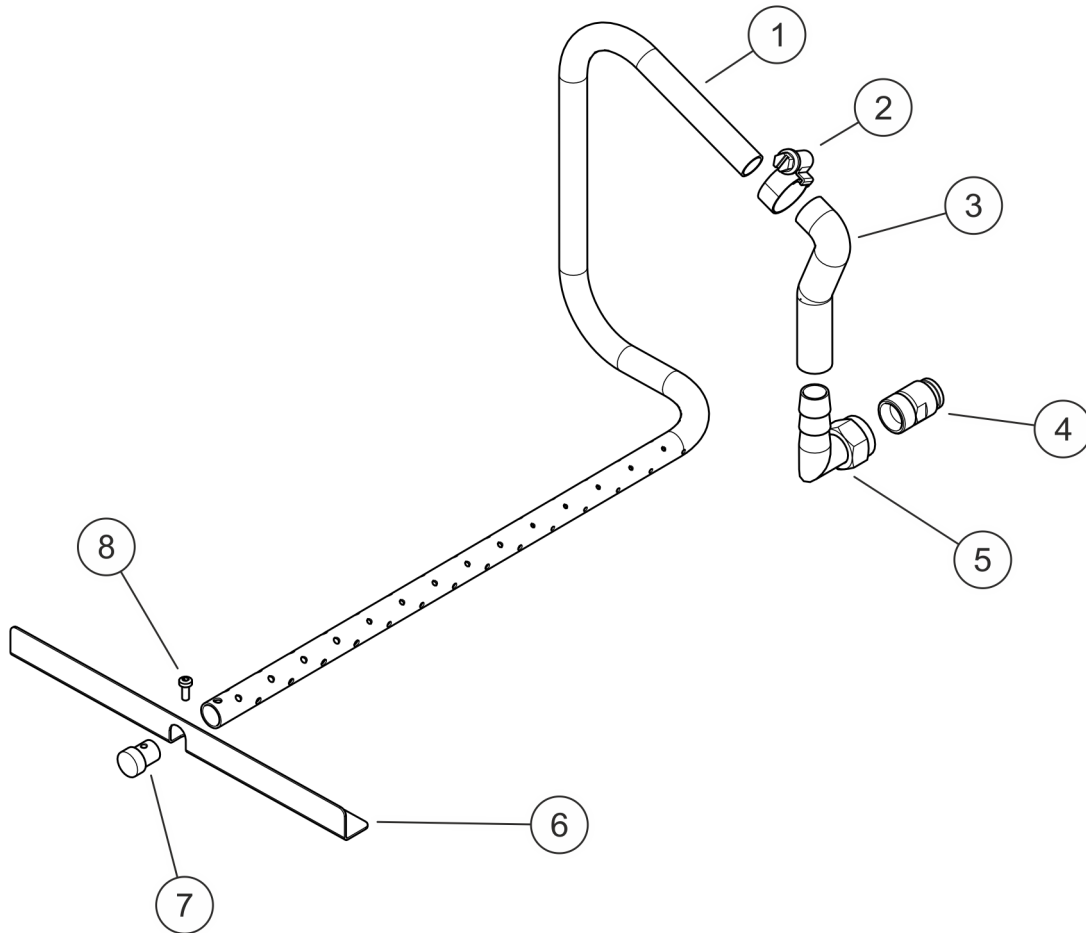
CAUTION!
Contact with hot / cold device parts, accessories and heat transfer liquid

Scalding, hot or cold burns

- Allow device parts, accessories and heat transfer liquid to reach room temperature before touching.

3 Scope of delivery

No.	Quantity	Description
1	1	Jet pipe
2	1	Hose clip
3	1	Shaped hose piece
4	1	Threaded nipple
5	1	Angle connector
6	1	Brace for jet pipe
7	1	Plug
8	1	Flat headed screw M3 x 08 (TX10)
-	1	Mounting instructions


4 Preparations


- | | |
|-----------------------|-----------------------------------|
| Personnel: | ■ Specialized personnel |
| Protective equipment: | ■ Protective work clothing |
| | ■ Safety gloves |
| | ■ Safety shoes |
| Tool: | ■ Allen key 4 mm |
| | ■ Crosshead screwdriver PZ2 |
| | ■ Crosshead screwdriver PH2 |
| | ■ Hexagon socket screwdriver TX10 |
| | ■ Open-end wrench 14 mm |
| | ■ Open-end wrench 19 mm |
| | ■ Water pump pliers |
| | ■ Socket wrench screwdriver 7 mm |

Mounting notes

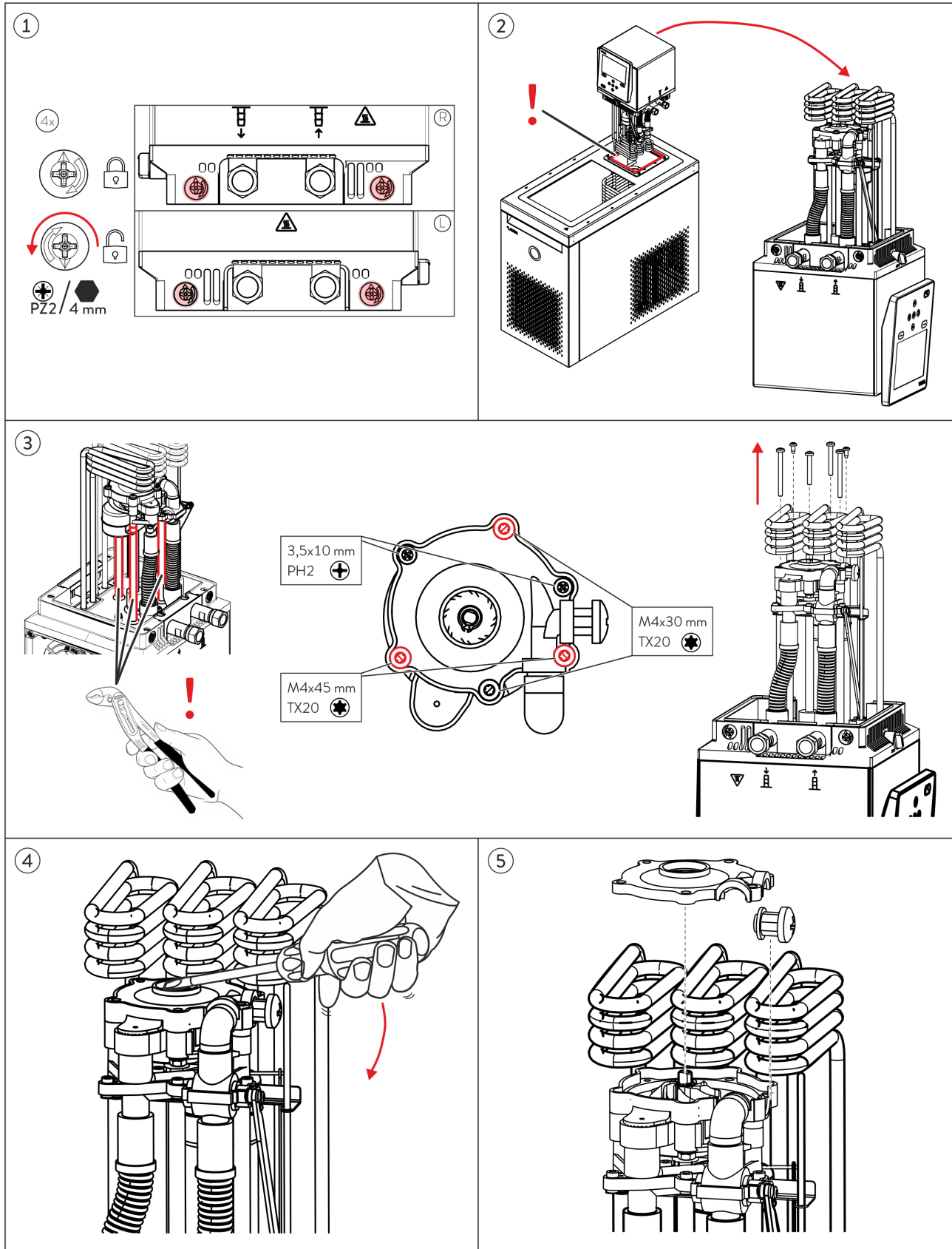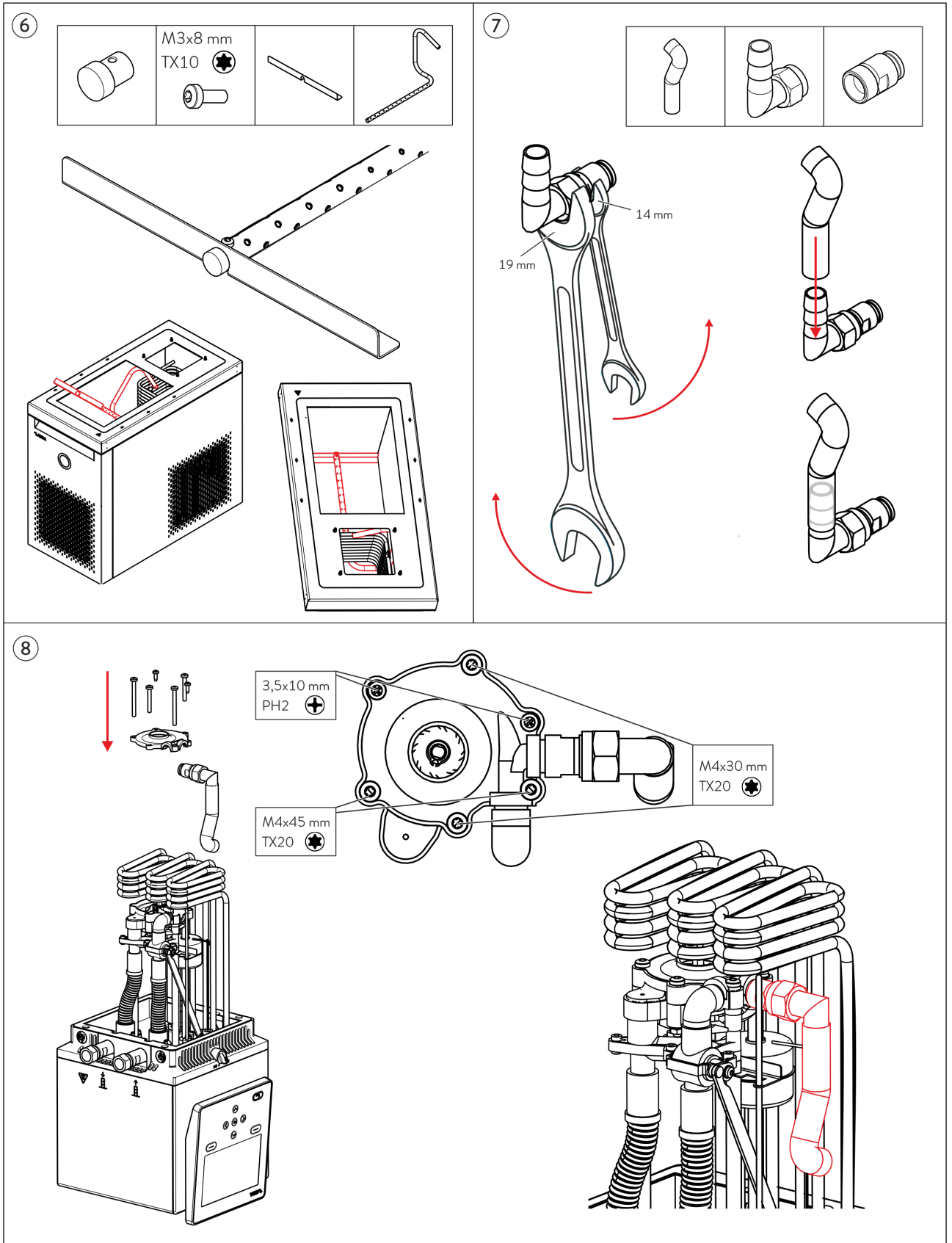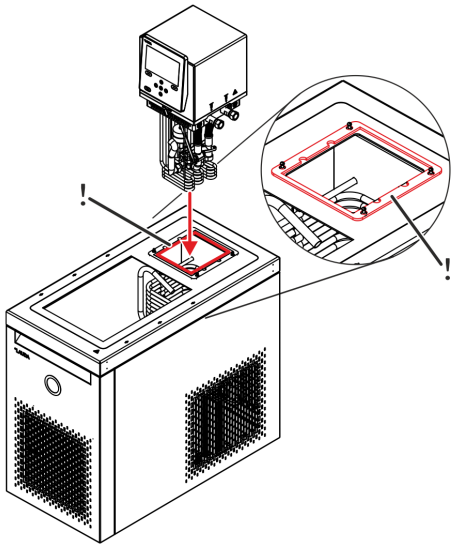
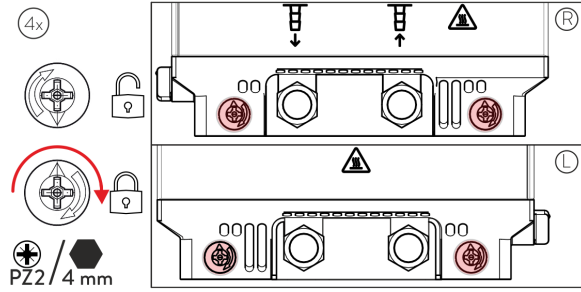
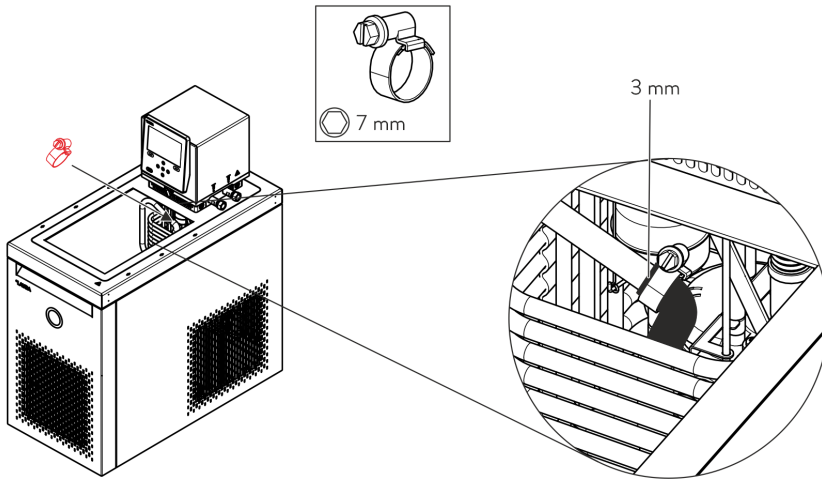
- Ensure that the device is disconnected from the power supply and secured against being switched back on.
- Ensure that the thermostat bath is completely empty and that the bath and the pump and control unit are clean and dry.
- Place the device on a stable surface and ensure there is sufficient space to place the dismantled pump and control unit.
- Check the scope of delivery of the accessory set for completeness and possible damage.
- No additional sealing material is required for the screw connections. The screw connections are tight if the contact surfaces are clean and undamaged and if they are assembled correctly.
- The shaped hose piece is secured to the jet pipe with one hose clip supplied with the accessory set. It is only pushed onto the angle connector.
- An FKM flat gasket is inserted on the bath bridge in the area of the locking bolts. When attaching the pump and control unit, ensure that the gasket is correctly seated between the bath bridge and the pump and control unit.
- **Please note the following regarding point ③ of the mounting instructions:**
Three screws are inserted from the underside of the pump chamber, through the pump chamber, and screwed into spacer columns above the pump chamber. When loosening these screws on the underside of the pump chamber, the spacer columns above the pump chamber must be held with a water pump pliers and secured against rotation.

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Q4DT-P_09-108, 1, DE_EN 09.04.2026 © 2025